


Control Standard Endotoxin

Instruction for use

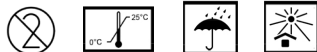

REF: ET-120V

Product Name: Control Standard Endotoxin

Specifications:

| Item No. | Product Name | Spec. |
|----------|----------------------------|---|
| ET-120V | Control Standard Endotoxin | E. Coli strain : O111: B4, 51-200 EU, 10 pcs/box |

Intended use:

Recommended for recheck the sensitivity of TAL, interference tests, and positive control.
It is intended for use as the control standard endotoxin in gel-clot and kinetic turbidimetric assay.

Notes:

Reconstitute the Endotoxin standard with BET water, the volume please refer to CoA accordingly,
Vortex for 15mins, for each dilution, Vortex 15s.

After reconstitution before any dilution, the CSE can be sealed and stored in refrigerator at -10C for 15days. When you need to use it, take it out and unfreeze it at room temperature and then Vortex for 15min.



Zhanjiang Bokang Marine Biological co., Ltd.

#188, RuiYunNan Rd., MaZhang district, Zhan Jiang, Guang Dong, China

Zipcode: 524094